LED-Kaltlichtquelle  
LED Cold Light Source

# LED3 Series

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Vielen Dank

# General Information

This high quality product has been designed and developed to require a minimum of maintenance. Please follow the guidelines below and read this manual in its entirety to ensure safe operation of the unit.

## Symbols

The following symbols are used in this manual.

CAUTION! This symbol indicates a potential risk and alerts you to proceed with caution.

## Warning

- Please always disconnect mains plug before cleaning the device.
- For indoor use only.

## Safety Information

- To avoid eye damage, don't look directly into the light source when turned on

**Important:** The evaluation of eye safety occurs according to the standard CIE S009/E:2002 („photo biological safety of lamps and lamp systems“). Within the risk grouping system of this CIE standard, the LEDs specified in this data sheet fall into the „low risk“ group (relaying to devices in the visible spectrum with an exposure time of 100s). Under real circumstances (for exposure time, eye pupils, observation distance), it can be assumed that no endangerment to the eye exists from these devices. As a matter of principle, however, it should be mentioned that intense light sources have a high secondary exposure potential due to their blinding effect. As is also true when viewing other bright light sources (e.g. headlights), temporary reduction in visual acuity and afterimages can occur, leading to irritation, annoyance, visual impairment, and even accidents, depending on the situation.

## Intended Use

Under the motto „ more light, less costs “ the LED3 can be perfectly used for most technical applications such as quality assurance, laboratories, industrial endoscopy (machine vision) and a lot more. Furthermore the LED light source is a cost efficient and energy-saving alternative to conventional cold light sources. The LED3 transfers a much larger part of the energy to visible light and it has a life span of roughly 50.000 working hours, consuming only about 36W.

The LED3 can run through the whole period in continuous operation, without additional costs and downtimes. The device is almost maintenance-free. The light source impresses by a modern and state-of-the-art design which also stands out by its modularity and its flexibility. The light source is passively cooled and thus amongst others perfectly suitable for cleanrooms. It can be positioned in a horizontal and vertical way.

# Function and Operation

## Package Contents

Cold Light Source LED3	1
Power supply unit	1
User Manual	1

## Technical Specifications

Weight	approx. 3,7 kg
Dimensions	259 × 174 × 81 mm (L × W × H)
Operating voltage	15 V DC, protection class III (safety extra-low voltage SELV)
LED life span	50.000 h (according to manufacturer)
LED	30 W power LED
Power consumption	approx. 36 W
Light guide access	standard Ø 15 mm
Active diameter of the light guide	max. 9 mm
Light control	analogous brightness control
Material	aluminium anodized
Colour	titan, grey and orange
Operating temperature	10°C to 35°C
Relative humidity	max. 95%, non condensated
Cooling system	passive cooling without fan

## Models

	LED3-P 100-010055	LED3-PC 100-010056	LED3-N 100-010057	LED3-W 100-010058
Luminous flux (light guide 1 m × Ø 9 mm)	ca. 730 lm	ca. 590 lm	ca. 630 lm	ca. 630 lm
Color temperature	5.700 K	5.700 K	4.000 K	3.000 K
Colour rendering index (CRI)	70	90	80	80

## Getting Started

The LED3 is a plug & play device very easy to use.


Please check before use whether...

...the mains plug is connected.


...the device is switched on (the blue LED light glows).

...the brightness is set to the desired value.

## Description of Buttons and Switches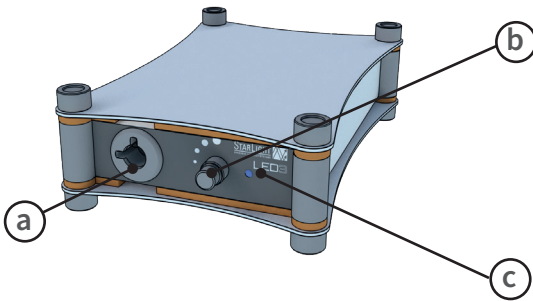

- a) **Light Guide Access**  
Connection of the light guide with a standard diameter of 15mm
- b) **Brightness Control**  
For the adjustment of the light (standby until 100%)
- c) **Operation Indicator**  
Blue LED lights when the device is turned on
- d) **Connection to Power Adapter**

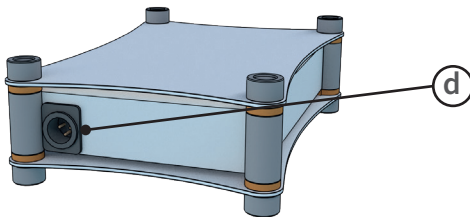# Technical Assistance

## Troubleshooting

The information in the table below can help to resolve the possible operating problems quickly.

Problem	Causes	Solution
No function at the device No blue LED	no network connection	Please check whether the power adaptor is connected.
	electronic defective	Please contact your distributor.
No light, blue LED glows	brightness adjustment on 0	Please adapt light to the desired value.
	LED and / or electronic defective	Please contact your distributor.
LED gets dark	thermal overload of the system	Please ensure an ambient temperature from 10° to 35°C.
	LED life span exceeded	Please contact your distributor.

## Storage, Maintenance, Repair

The device is maintenance-free. Some points, however, you should take into consideration.

1. Please unplug the device before cleaning.
2. For the cleaning of the housing please use a smooth, semi-wet cloth, if necessary with a smooth household cleaning fluid.
3. No moisture should entering into the housing interior.
4. To avoid damages during storage or transport please retain the packaging.

**Repairs are only allowed by trained and experienced personnel. In case of improper handling the warranty expires.**

## Technical Service

### Web Resources

Visit our StarLight website at [www.starlight-sl.de](http://www.starlight-sl.de) for:

- Contact details of technical customer service.
- StarLight product overview and price list as well as infos about accessories and spare parts.
- Additional product information and special offers.

### Contact

For information or technical assistance contact your local distributor or directly StarLight Sensorik - Lichtleitsysteme GmbH & Co. KG.

## Warranty

StarLight Sensorik - Lichtleitsysteme GmbH & Co. KG warrants that this product will be free from defects in material and workmanship for a period of one (1) year from date of purchase. If a defect is present, StarLight will, at its option, repair, replace, or refund the purchase price of this product at no charge to you. Precondition is the device is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, or misapplication, or from ordinary wear and tear.

For your protection, items being returned must be insured against possible damage or loss. This warranty shall be limited to the replacement of defective products.

It is expressly agreed that this warranty will be in lieu of all warranties of fitness and in lieu of the warranty of merchantability.

## Equipment Disposal

This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste. Instead, it's your responsibility to correctly dispose of your equipment at lifecycle -end by handling it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

Thank you

## Notizen / Notes